


PAINTED STEEL WORK SURFACES

PRODUCT DEFINITION

Teclab painted steel tops are made from furniture grade cold rolled steel. Features include welded corners, reinforcing channel members that stiffen the entire work-surface and a durable powder coated finish to match or compliment your casework.

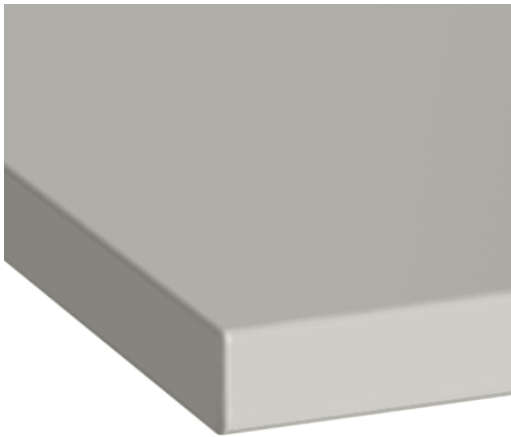

PRODUCT SPECIFICATIONS


Thickness - 1"

Material - 18-gauge cold rolled steel

Finish - Textured powder coat paint

Sizes - Up to 46" x 96"

Edge - Wrapped and welded corners

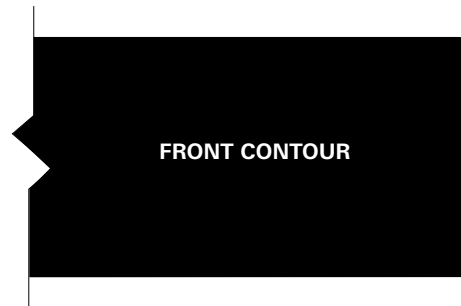

GENERAL CARE & MAINTENANCE

Powder coated steel is a chemical resistant, baked coating that is able to withstand a number of chemicals. However, in order to prolong the "like new" appearance of your casework installation, it is necessary to regularly clean all exposed surfaces.

Regular cleaning

Powder coated steel casework needs to be cleaned regularly to remove surface contaminants. This also helps to safeguard against unwanted chemical reactions. The best method for cleaning is to use glass cleaner with a soft cloth. Be sure that this cleaner is compatible with other chemicals that may be in use around the area you are cleaning as ammonia based products should never be mixed with products containing bleach.

Caution:


- **Do not** use steel wool or abrasive cleaners for cleaning. This may damage the surface(s) of your cabinet(s).
- **Do not** expose powder coated surfaces to acids which may cause staining or damage to the exposed surface(s).

Chemicals to Avoid:

- EK (Methyl Ethyl Keytone), Acetone, or other strong solvents
- Chlorine-based cleaners
- Sulfur-based cleaners

FINISH OPTIONS

Powder coated steel.

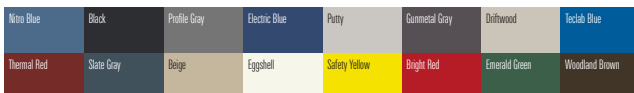

Custom paint color matching is available.